





Each punch tip is manufactured to a very high degree of accuracy. Even a slight wear & tear can result in production of dull or collared tablet. This makes polishing the punch cavity at regular interval an important process to ensure consistent tooling performance.

The Punch Cavity Polishing Kit helps you maintain the original cavity finish. This kit contains different types of brushes for different cavities & different types of pastes for varied application.

Polishing is done to revive the original cavity finish & not to remove material. Hence precaution needs to be taken to prevent excessive polishing. The 4 most common errors in punch cavity polishing are:

- a. Rounding of the outside diameter of the tip
- b. Rounding of the edges of the tip face
- c. Distortion of embossed details
- d. Distortion of the tip face

KIT CONTENTS

- · Polishing Motor with 1/4" Chuck
- Felt Bob Dome Type
- Felt Bob Conical Flat Type
- . Nylon Brush Cup Type
- Nylon Brush Wheel Type
- · Diamond Paste
- . Cavity Shining Paste
- Hiffin Lubricant
- Emery Flour Powder

TECHNICAL SPECIFICATIONS


Motor : 1/8 HP, 2800 RPM

Electrical Supply : 110V/1PH OR 220V/1PH

Dimensions : 50 x 50 x 25cms

Net Weight : 10 kgs

Gross Weight : 14 kgs

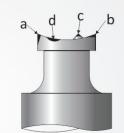







*specifications subject to change without notice



52, Suyog Industrial Estate, LBS Marg, Vikhroli (W), Mumbai - 400 083, India. T: +91(22) 25786007, 25793111 F: 25771900 E: pi@pharmachine.com W: www.pharmachine.com



www.facebook.com/TabletPressAdeptEngineers www.adepttabletpress.blogspot.in www.slideshare.net/AdeptTabletPress

www.youtube.com/user/AdeptTabletPress